

How to activate simple software products on Software Portal

The following products may be activated on <https://activation.mt.com>


1. LabX
2. Calibry
3. MC Link
4. ProdX
5. Collect+
6. Form+
7. EasyDirect PAM, from 2.0
8. PipetteX

→ upload License key file **A**

9. EasyDirect pH, from V1.1
10. EasyDirect DE/RE
11. EasyDirect Titration, from V1.3
12. EasyDirectMoisture
13. QuickPredict Embedded Moisture
14. EasyDirect Balance
15. Data+
16. LabClimate

→ enter License key **B**

If you have type **A** license keys, please refer to the link "[Activation Manual E-Learning](#)" on the Main screen for instructions:

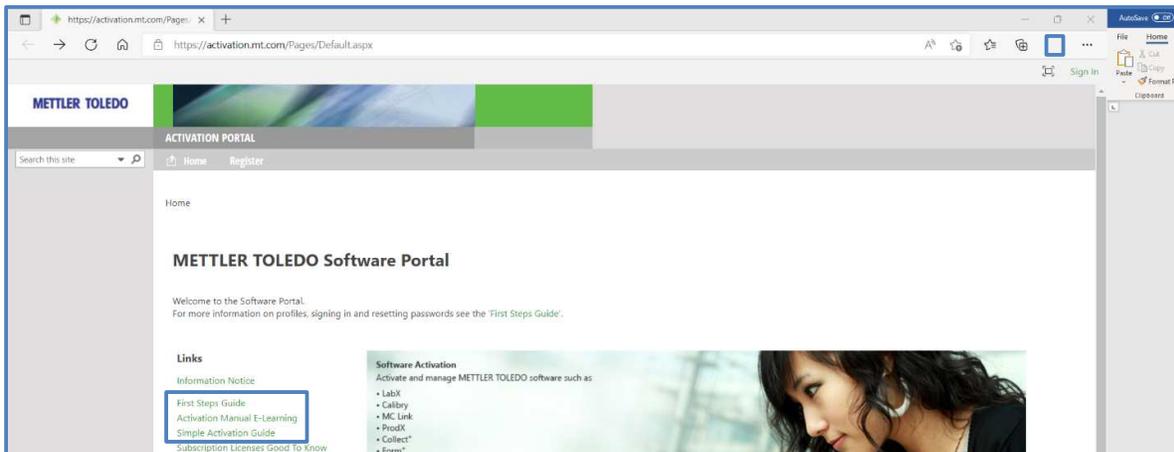


This guide describes how to activate a type **B** license key. Preconditions:


- You have an installed type **B** SW product.
- You have a license key for the same type **B** product.
- You have registered for a user account. Please refer to the [First Step Guide](#) on the main screen for instructions.

Steps to activate a type **B product:**


1. Start your installed type **B** software product and note / copy the License Registration Information (LRI)

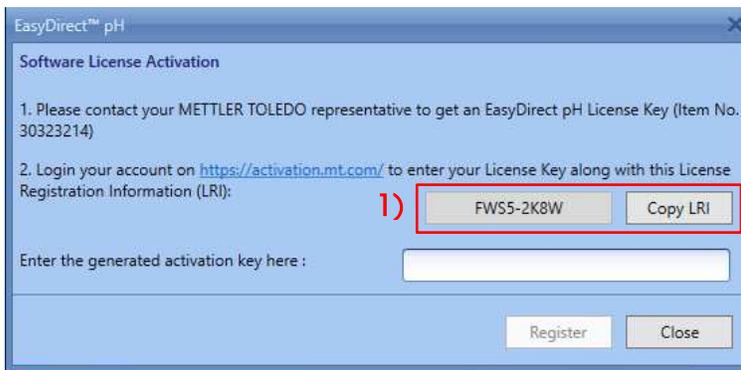

2. Sign in to <https://activation.mt.com> and refer to "Activation"
3. Enter your type **B** license key and enter / paste the LRI and submit

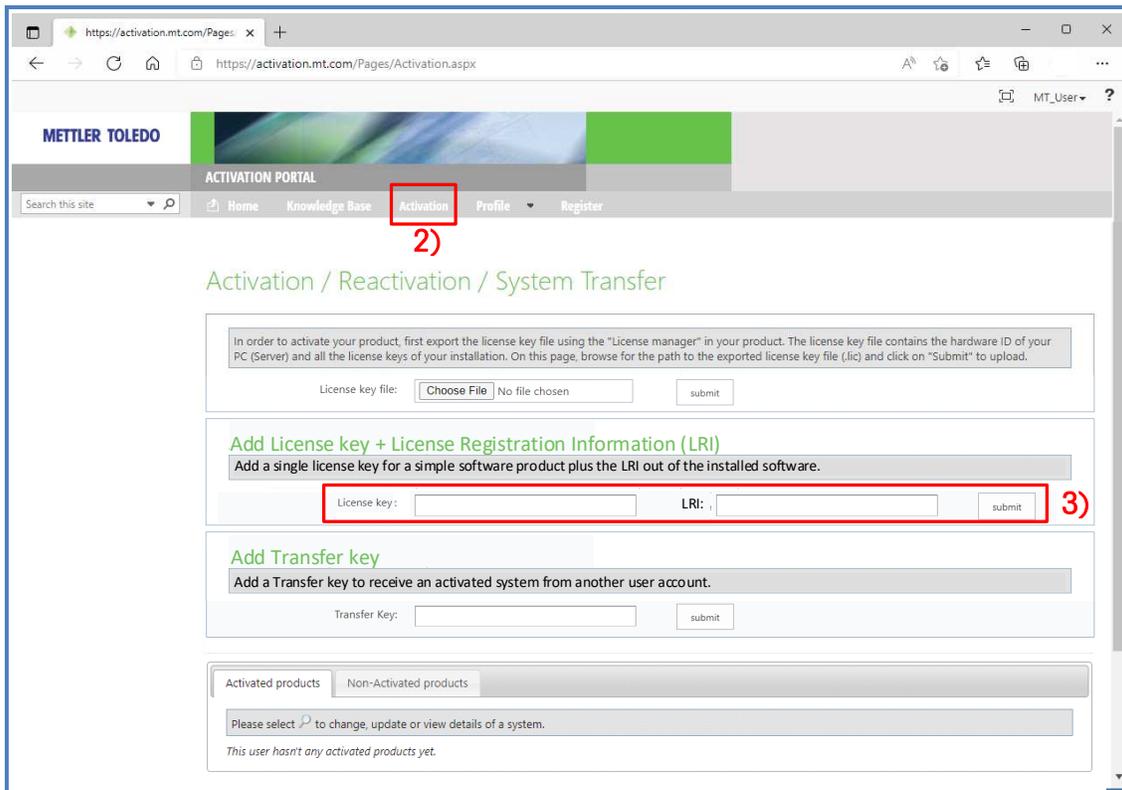

- The Software Portal provides a Short Activation Code (SAC) in a pop-up box.
Note / copy the Short Activation Code (SAC)

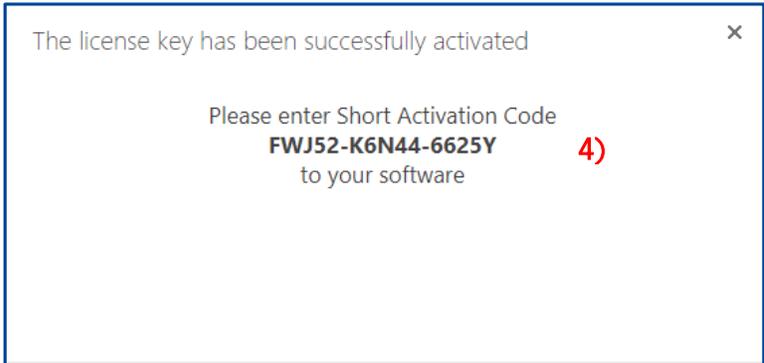

- Enter / paste the Short Activation key (SAC) into your Type **B** software and register.



Your type **B SW is now activated!**

View your activated system on Software Portal:

The screenshot shows the Mettler Toledo Activation Portal interface. The browser address bar displays <https://activation.mt.com/Pages/Activation.aspx>. The navigation menu includes Home, Knowledge Base, Activation (highlighted with a red box), Profile, and Register. The main heading is "Activation / Reactivation / System Transfer".

Below the heading, there are three main sections for activation:


- License key file:** A file upload section with a "Choose File" button and a "submit" button.
- Add License key + License Registration Information (LRI):** A section for adding a single license key and LRI, with input fields for "License key:" and "LRI:" and a "submit" button.
- Add Transfer key:** A section for adding a transfer key, with a "Transfer Key:" input field and a "submit" button.

At the bottom, there is a table titled "Activated products" (highlighted with a red box). The table has two tabs: "Activated products" and "Non-Activated products". The table contains one row of data:


Product Subname	System name	Server name	Server location	Hardware ID / LRI
EasyDirect pH				FW55-2K8W

Add / edit system information (optional)

1. Select an activated system and click the magnifier

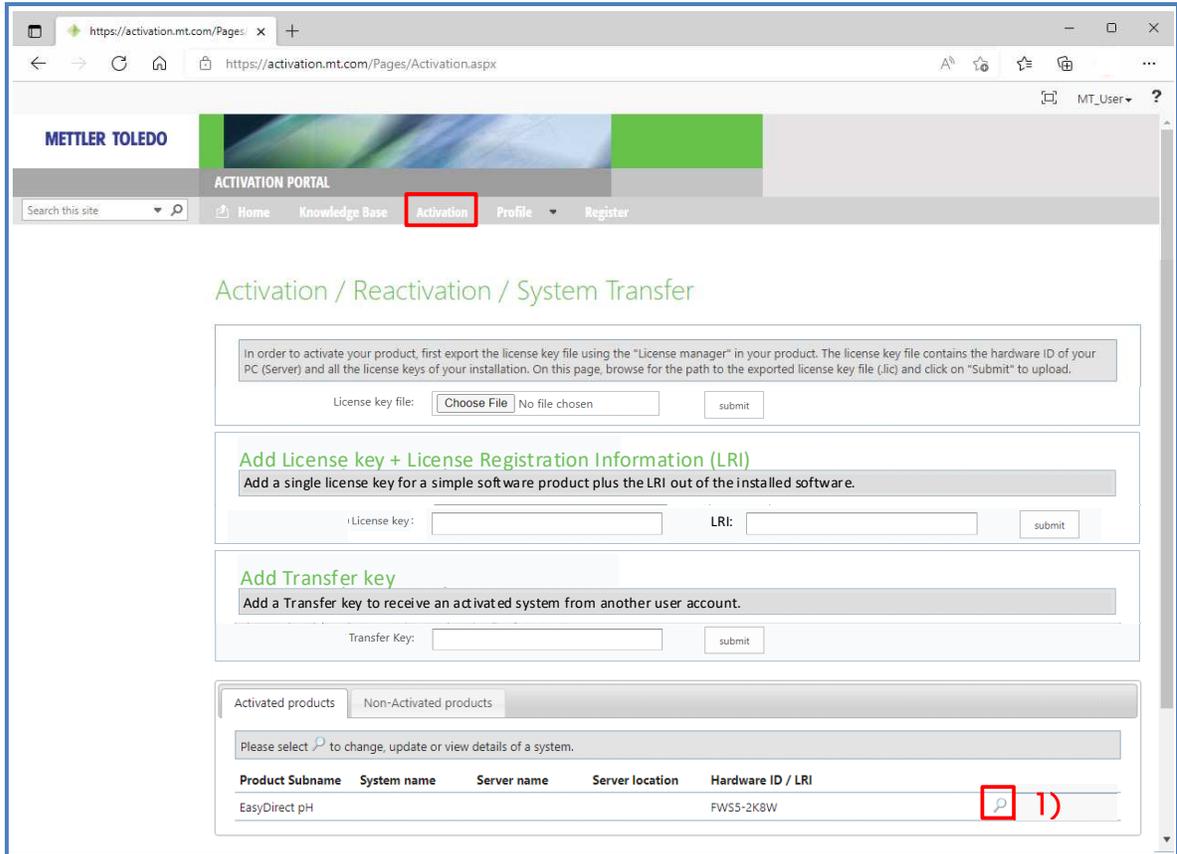

2. Select Edit System in License overview and add/change system information

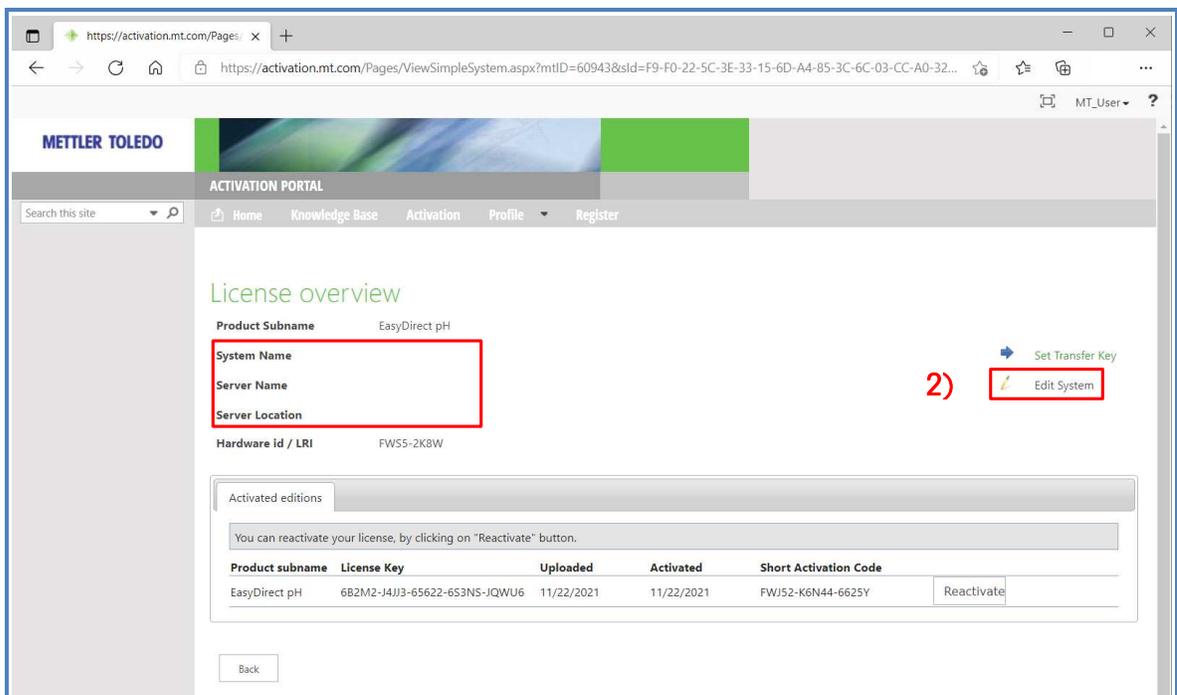

3. View activated systems including system information

The screenshot shows the Mettler Toledo Activation Portal interface. The navigation menu includes 'Home', 'Knowledge Base', 'Activation', 'Profile', and 'Register'. The 'Activation' menu item is highlighted with a red box. The main content area is titled 'Activation / Reactivation / System Transfer' and contains several sections for license management: 'License key file' with a 'Choose File' button, 'Add License key + License Registration Information (LRI)', and 'Add Transfer key'. At the bottom, there is a table of 'Activated products' with a search icon and a red circled '3)' next to it. The table has columns for Product Subname, System name, Server name, Server location, and Hardware ID / LRI. The first row shows 'EasyDirect pH' with system name 'pH-31', server name 'PCC 337', server location 'Room 4-203', and hardware ID 'FW55-2K8W'. The 'System name', 'Server name', and 'Server location' columns are highlighted with a red box.

Product Subname	System name	Server name	Server location	Hardware ID / LRI
EasyDirect pH	pH-31	PCC 337	Room 4-203	FW55-2K8W