

How to activate simple software products on Software Portal


The following products may be activated on <https://activation.mt.com>

1. LabX
2. Calibry
3. MC Link
4. ProdX
5. Collect+
6. Form+
7. EasyDirect PAM, from 2.0


→ upload License key file **A**

8. EasyDirect pH, from V1.1
9. EasyDirect DE/RE

→ enter License key **B**

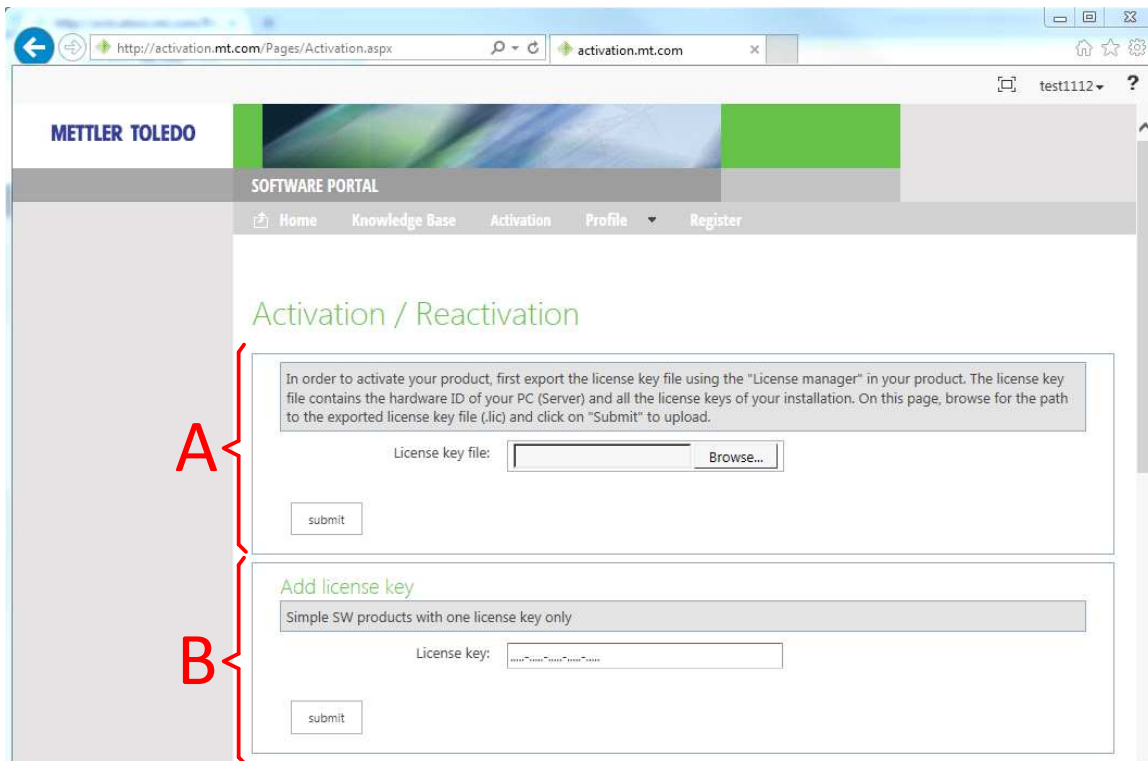

If you have type **A** license keys, please refer to the link "Activation Manual E-Learning" in Main screen for instructions:

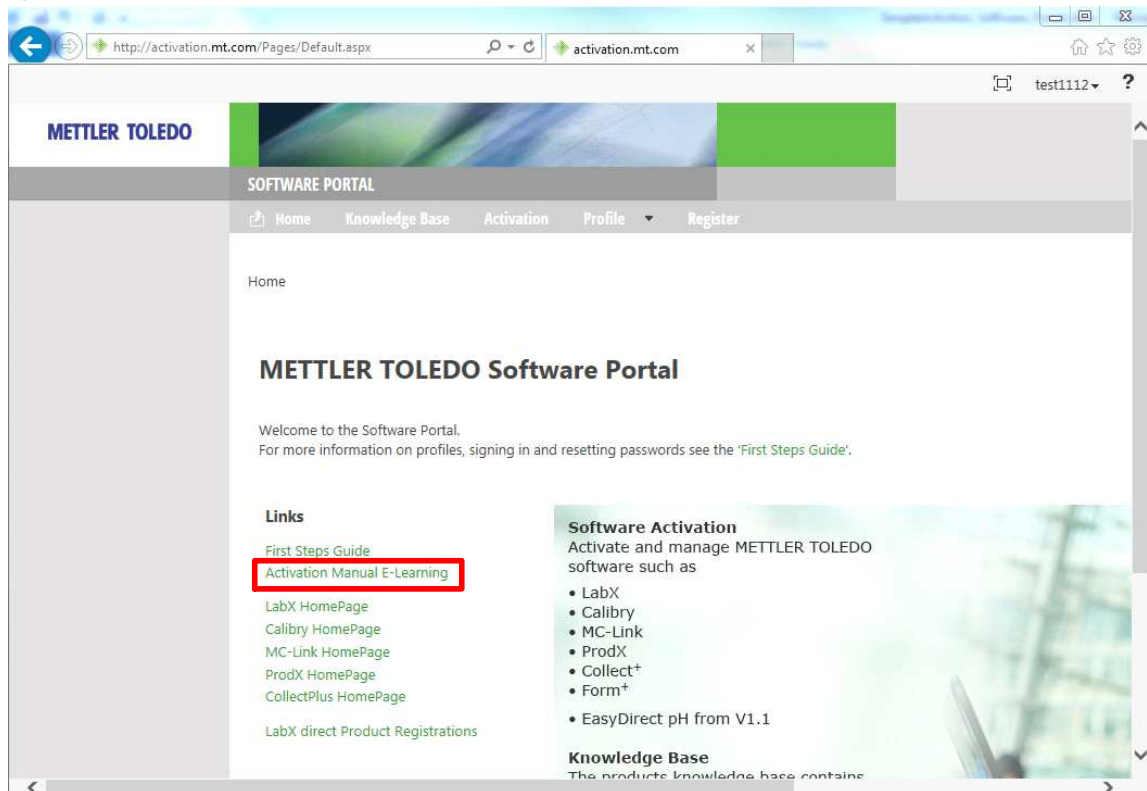


This guide describes how to activate a type **B** license key.

Preconditions:


- You have a type **B** SW product installed on a PC
- You have a license key from a type **B** product

Steps to activate a type **B** product:


1. Sign in to Software Portal
2. Refer to "Activation"
3. Enter and submit your type **B** license key

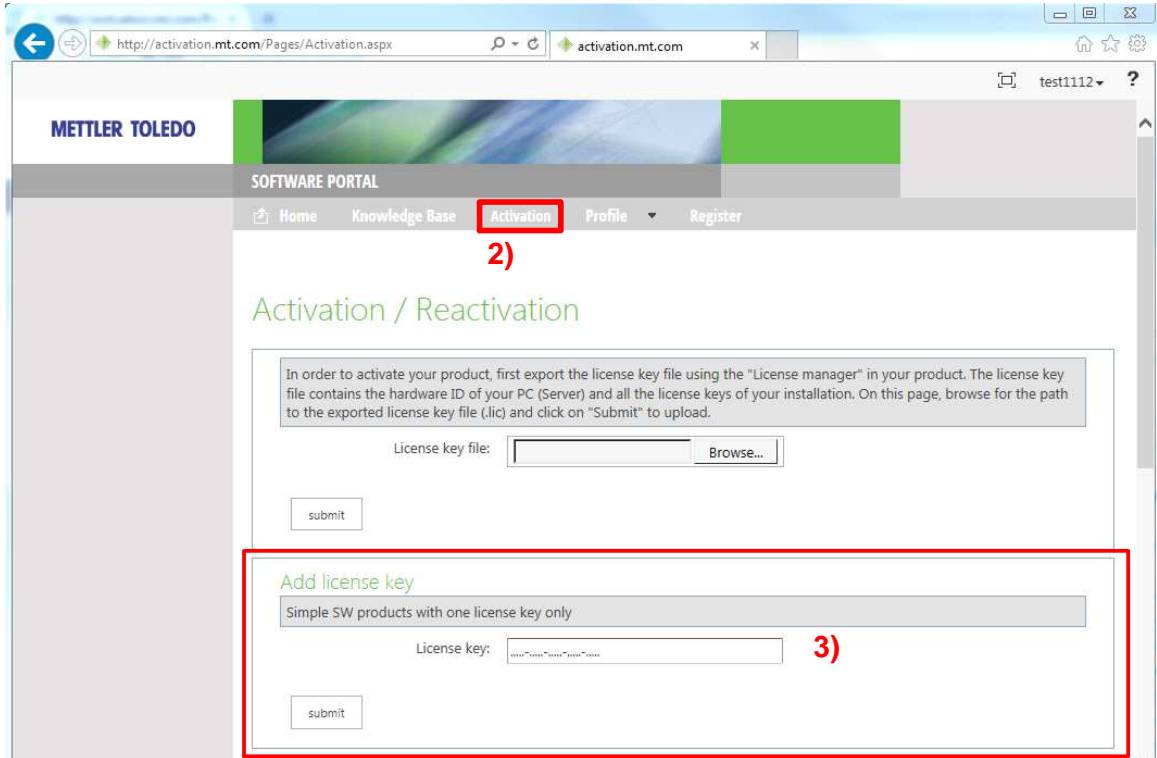


4. Refer to the type **B** software product, installed on the PC

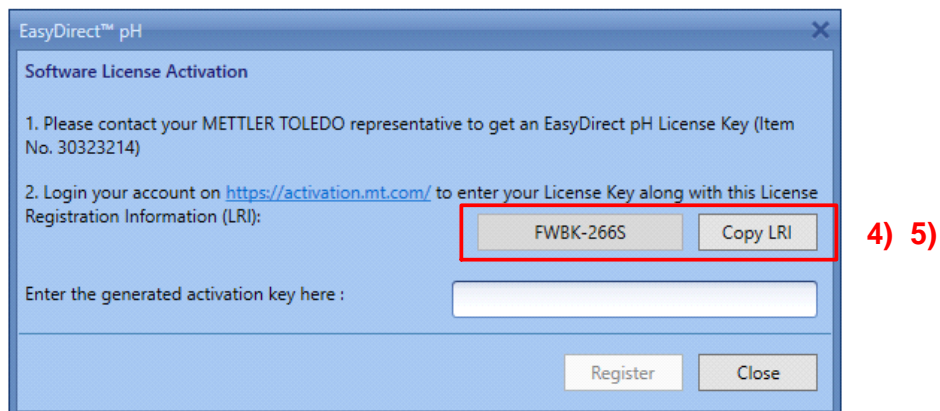


- 5. Copy License Registration Information (LRI)
- 6. Add the LRI to the previously uploaded license key and

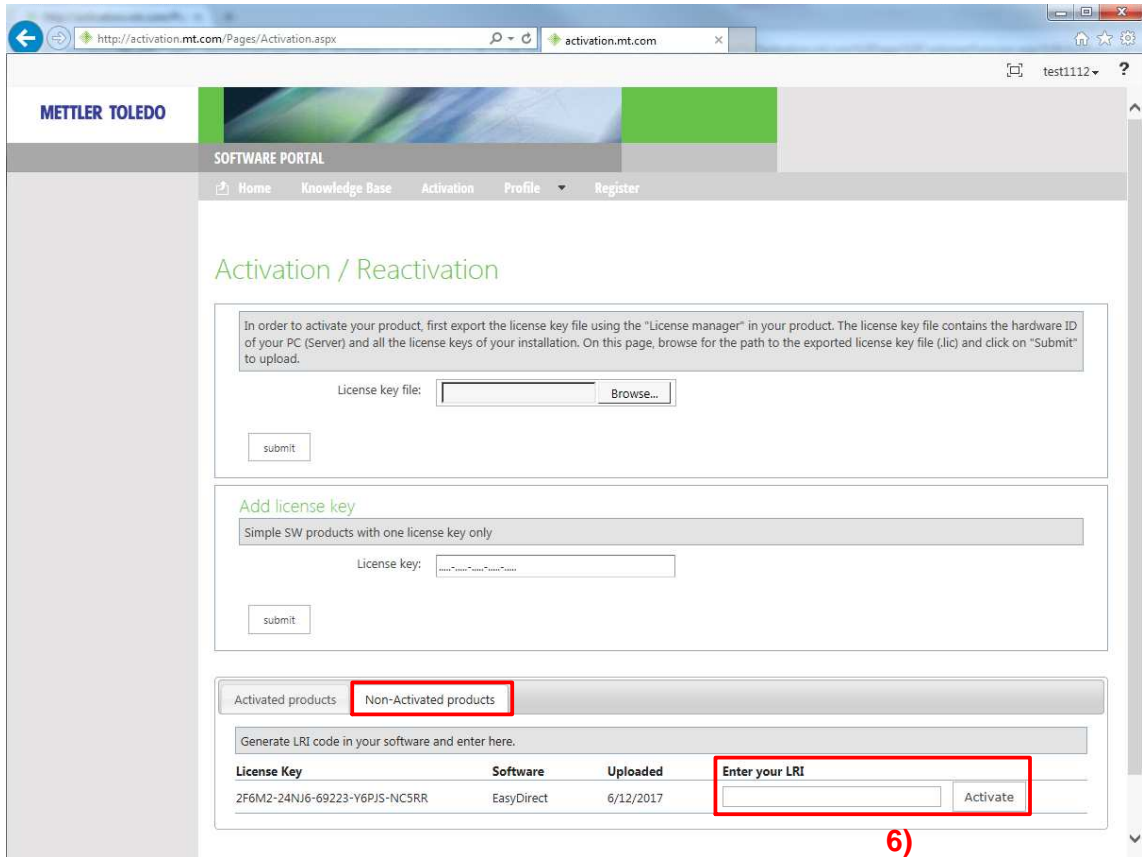


- 7. Add system information, at least a system name

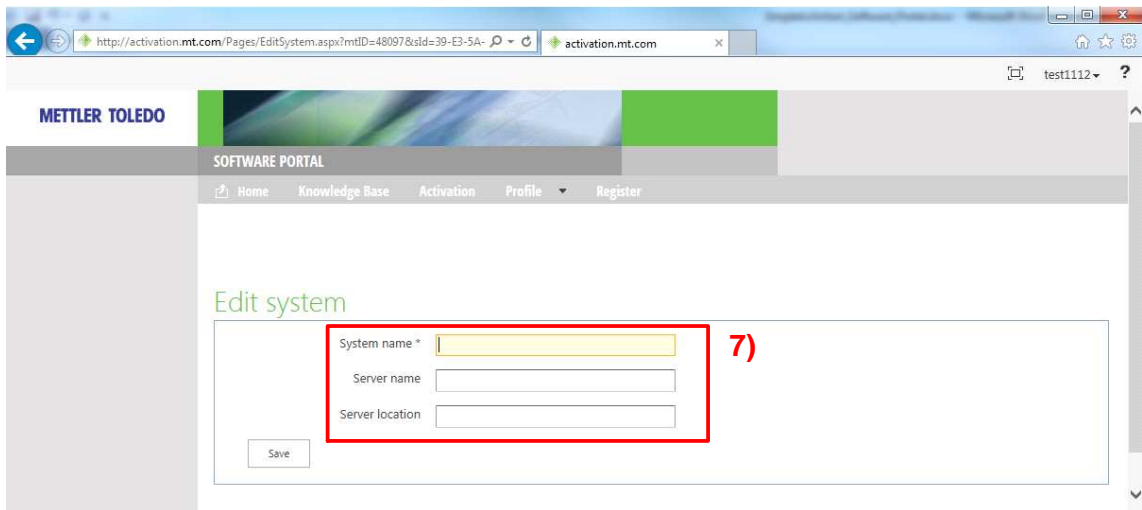

8. Copy Short Activation Code (SAC) from pop-up box

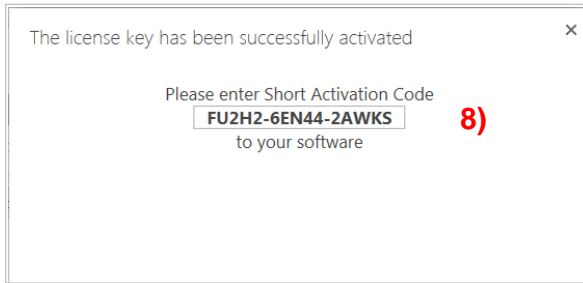

9. Enter the Short Activation Code into the type **B** SW installed on PC

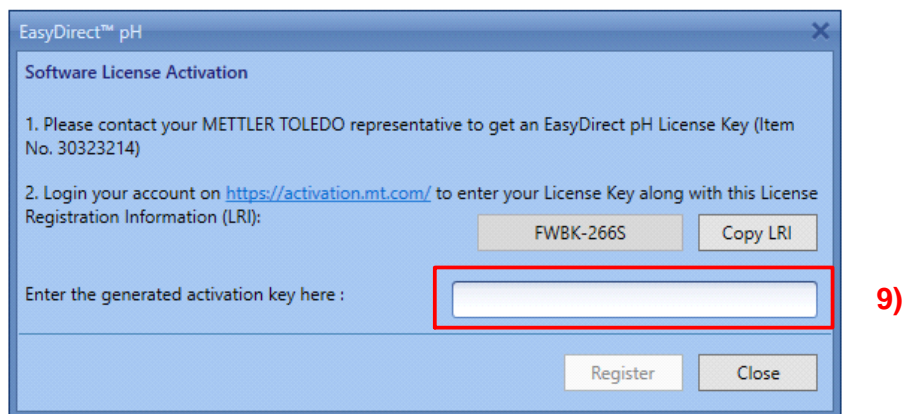

10. Your type **B** SW is now activated!

11. View your activated systems on Software Portal

The screenshot shows the 'Activation / Reactivation' page on the Mettler Toledo Software Portal. The page includes a navigation menu with 'Home', 'Knowledge Base', 'Activation', 'Profile', and 'Register'. The 'Activation' tab is selected. Below the navigation, there are two main sections: 'License key file' and 'Add license key'. The 'License key file' section has a text box and a 'Browse...' button. The 'Add license key' section has a text box and a 'submit' button. Below these sections is a table of activated products. The table has columns for 'Product name', 'System name', 'Server name', 'Server location', and 'Hardware ID / LRI'. The first row is highlighted with a red box and the number 11.

Product name	System name	Server name	Server location	Hardware ID / LRI
EasyDirect	Test201706121337			FSAD-266N
LabX	Test201706111144			ALL6G-VKF2L- YE63F-F6S8K-ERFRX
EasyDirect	Test201706111138			BYKS-266N
EasyDirect	Test201706111100			FSAD-266N
LabX	Test201705041437			ALL6G-VKF2L- YE63F-F6S8K-ERFRV