

Subscription licenses

good to know

SW products with subscription licensing

The following product(s) will expire after a certain period:

1. EasyDirect PAM from V2.0

License types:

Edition licenses

- Starter
- Site
- Unlimited

Subscription Edition licenses unlock a certain number of functions, depending on the Edition level (Starter, Site, Unlimited) for a duration of 12 month.

Renewal licenses

- Starter renewal
- Site renewal
- Unlimited renewal


Subscription Renewal licenses extend the according Edition license functionality for a duration of 12 month.


A renewal license can only get activated, if a valid license of the same level has been activated previously. Valid pre-activations are:

- For Starter renewal: Starter edition or Starter renewal
- For Site renewal: Basic edition or Basic renewal
- For Unlimited renewal: Unlimited edition or Unlimited renewal

Expiry date

As soon as a new license has been activated, the expiry date is shown in the System:

- 1. Select Activation tab in menu
- 2. Select system 
- 3. View expiry date

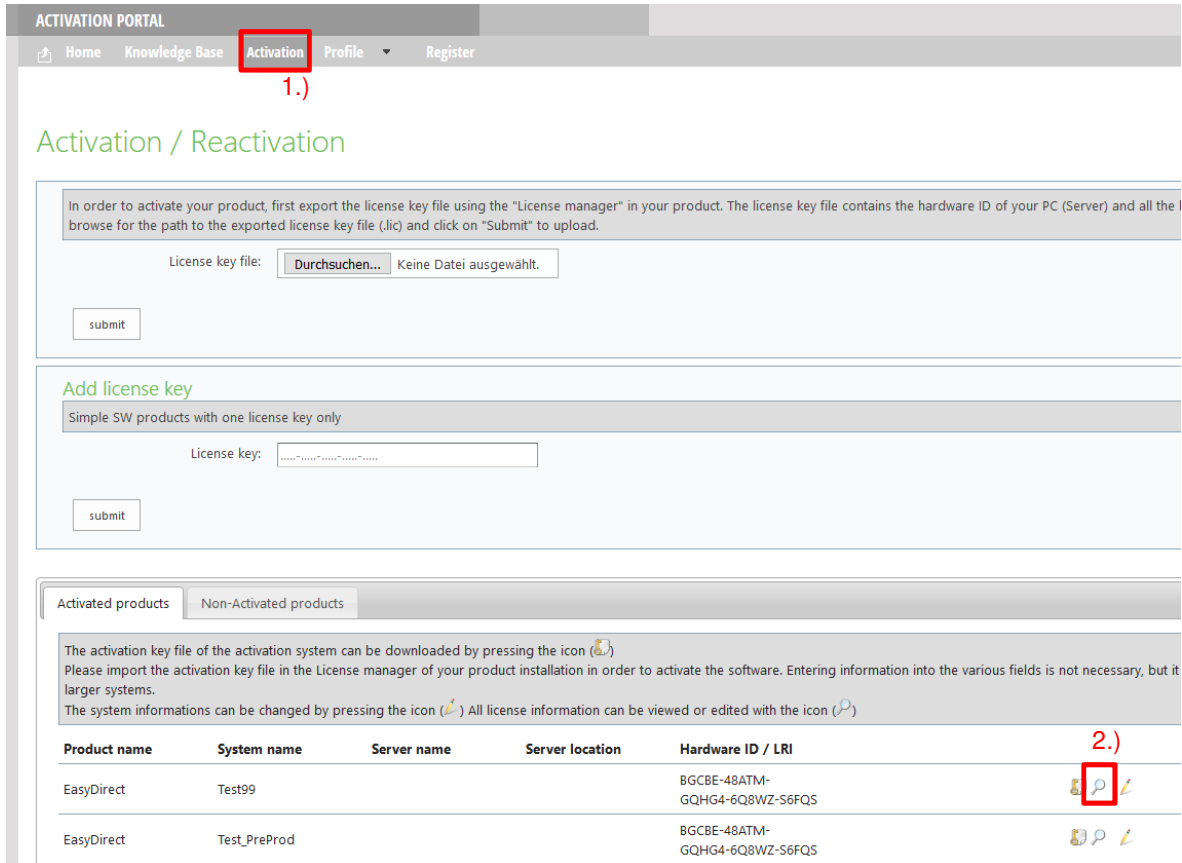

ACTIVATION PORTAL

Home Knowledge Base **Activation** Profile Register

1.)

Activation / Reactivation

In order to activate your product, first export the license key file using the "License manager" in your product. The license key file contains the hardware ID of your PC (Server) and all the...
browse for the path to the exported license key file (.lic) and click on "Submit" to upload.

License key file: Keine Datei ausgewählt.



Add license key

Simple SW products with one license key only

License key:

Activated products Non-Activated products

The activation key file of the activation system can be downloaded by pressing the icon (📄)
Please import the activation key file in the License manager of your product installation in order to activate the software. Entering information into the various fields is not necessary, but it...
larger systems.
The system informations can be changed by pressing the icon (✎) All license information can be viewed or edited with the icon (🔍)

Product name	System name	Server name	Server location	Hardware ID / LRI	
EasyDirect	Test99			BGCBE-48ATM- GQHG4-6Q8WZ-56FQS	
EasyDirect	Test_PreProd			BGCBE-48ATM- GQHG4-6Q8WZ-56FQS	

2.)

ACTIVATION PORTAL

Home Knowledge Base Activation Profile Register

License overview

Edition and option licenses

Edition and option licenses for test 2345 (Test99)

Active licenses

Editions				
Title	License key	Type	Activation date	Expiry date
Pipette Asset Management 2.0 Enterprise	2H6S2-KJ6P6-27422-26KJS-2L727	sidegrade	4/5/2018	4/5/2019 3.)

Options
No registered license for this type.

Inactive licenses

Reminder email

60 days before a license key expires, <https://activation.mt.com> sends a reminder email to:

- The user email address
- A second email address, if a 2nd email address is stored in the user profile:

How to add a second email address:

- Select "Profile tab in menu"
- Select "View Profile"
- select "Edit"
- add the 2nd email address into "E-mail (secondary)"

As soon as you receive a reminder email, please contact your local METTLER TOLEDO representative, in order to order a renewal license. When the license is expired, the application will refuse to work.